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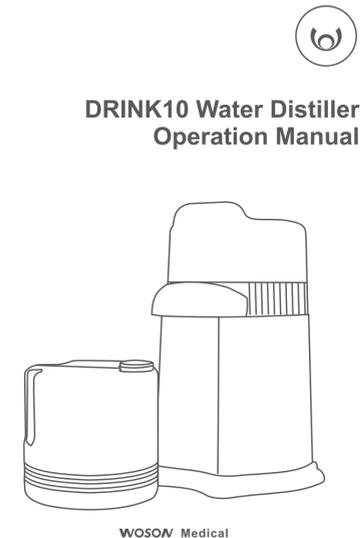
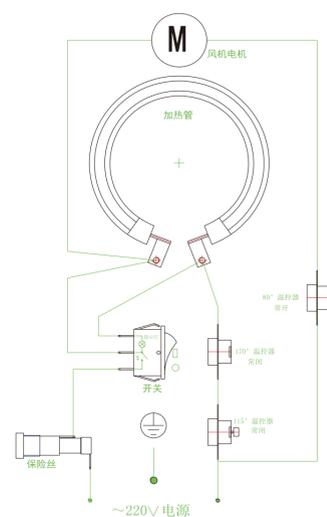
Chapter 3 Receiving and Installation

3.1 Checking the Package

Please check the package carefully when you receive the production, for the production can be damaged by package disrepair anywhere.

Part Name	Function
KSD301 Temperature Controller	170 degree C automatic reset, Protecting Electric heating rings from damage of high temperature.
KSD301 Temperature Controller	115 Degree C manual reset, Avoiding auto-heating after water boiled while power on.
KSD301 Temperature Controller	80 degree C automatic reset, For energy-saving and safety, the cooling fan's running is controlled by chamber temperature.

Part Name	Function
Fuse	Clear the circuit when short circuit or over current.
Rocker Switch (Power Switch)	Convenient to start or end the distiller at any time

CAUTION MANUFACTURER IS NOT HELD RESPONSIBLE FOR ANY ARBITRARY DISASSEMBLY, AMENDMENT WITH THE UNIT, BY UNAUTHORISED PERSON OR UNPROFESSIONAL TECHNICIAN

2.3 Safety Parts

For avoiding risk, please be cautious at operation.

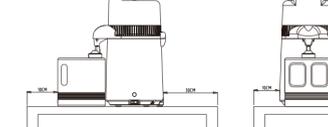
Operating Risk
For safety, please protect the cover after open it, for it can close automatically under external pressure.

Burn hazard
Do not open the cover while working, because you may be burned by water vapour.

The risk of contamination
For avoiding leaving contamination in chamber, please clean chamber as soon as possible.

3.3 Working Environment

Distiller installation should be left at least 10cm all around, and 20cm to the top.


Advise installing in an environment with good condition of ventilation.
Ambient temperature: 30℃.
Relative humidity: <80%
Air-pressure: 860hPa ~ 1060hPa
Earthing

3.4 Installation

- ✦ Set the distiller on a stable surface, and keep at least 10cm from graphic edge.
- ✦ For safety, please remove all stuff from distiller.
- ✦ For convenient operation, please remove all stuff in front of distiller.
- ✦ Do not set bottle with rusty liquid around distiller, for eliminating potential hazard while Thermal distillation.

3.5 Power Connector

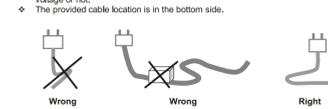- ✦ Check the nameplate under distiller bottom whether it corresponds relevant supply voltage or not.
- ✦ The provided cable location is in the bottom side.

Chapter 4 Product Description and Specification

4.1 Front View

No.	Name	No.	Name
1	Filter Tip	7	Water tank
2	List	8	Gasket ring
3	Body	9	Rubber Plug
4	Fuse Holder	10	Insulator Cover
5	Temperature Protection Switch	11	Chamber
6	Power Switch		

4.2 Size and Capacity

Size	DRINK10
Extensil Size	225*270*370mm
Packing Size	260*260*420mm
Full Level Capacity	Φ180*142

4.3 Specification

Essential Paramter
Rated Voltage: a.c.220V~240V/50/60Hz
Rated Power: 750VA
Fuse: 250v 1T10A
Operating Temperature: 5~40℃
Noise: < 50db
Speed: 1.2L/H
Water Rate: >= 96%

Chamber
Material: Stainless Steel304
Maximum Capacity: 4L
The Highest Temperature: 120 ℃
Volume Capacity Size: Φ180*185
Temperature Safety Protection:
Visual temperature protection: 115 ℃
Heating element temperature protection: 170 ℃

Chapter 5 Operation Process

5.1 Setting

Open the packing chest. Take the distiller out of the chest. Then take the water tank out of the distiller. Set the distiller on a stable surface in terms of picture 1.

5.2 Water-adding

Open the lid (As picture 2). Pour proper volume of water into the chamber within the highest waterline.
ATTENTION As the cover connects tightly with the body, please hold the body with one hand, and pull the cover upward with another.

5.3 Power Connector

Set the storage bucket stably according to picture 1, insert the one end of power cord into the body, and another into correct electric outlet.

5.4 Starting

Turn on the power switch (If the power switch off, press the heating button). If the indicator light on, it means start to heating. After a period of time, the cooling fan will run automatically for distilling.
It takes about 3.5 hours to distill 4L water, and the first pure drop will appear after half an hour, subject to local temperature.

5.5 Ending

When distilling finished, the heating system will stop automatically, and indicator light will be off. However, the cooling fan will keep running until the chamber temperature drops below 70 degree C.

CAUTION Heating button is in charge of heating process start. The heating process will stop automatically when there is no water or no water for distilling. (When heating, indicator light in power switch on; stopping, off). Power switch controls both of heating and cooling fan's running.
For filtering, the distiller is furnished with active carbon. In accordance with demand, you can put the active carbon into the faucet, and then pull it, you can take it off.

CAUTION As upper cover is heavy, do not open and erect it while using.

5.6 Power Off

Please turn off the power switch when no longer in use.
Please pull the end of power cord out of electric outlet when no longer in use.

Chapter 6 Maintenance

6.1 Maintenance Schedule Chart

Maintenance Required	Person Responsible
Daily	
Clean Body	User
Clean Chamber	User
Clean Water tank	User
One Month	
Clean Chamber with Abluent	User
One Quarter	
Filter Tip Replacement	User

6.2 Daily Maintenance

Please clean and drain the chamber and body with soft towel after using.
WARNING DO NOT CLEAN DOOR OR ASSEMBLE CHAMBER WITH METALLIC BRUSH, STEEL WOOL, OR PRODUCTS CONTAINING CHLORIDE. TO AVOID SCALD, PLEASE MAKE SURE DISTILLER IS COOLED DOWN BEFORE CLEANING.
Cleaning Water Tank
Water tank can be disinfected and cleaned with warm (not) water, and it will not be out of shape by hot water.

6.3 Monthly Maintenance

Cleaning Chamber with Abluent
The internal of stainless steel tank need to be washed with clean water in order to remove the distillation residual precipitation and impurities. And please clean the stainless steel tank per month (or depend on the quantity of dirt within the tank), with the detergent which we recommend to you. The procedure as follow:
A. Add enough (hot) water into tank till cover dirt.
B. Put one or four (depends on quantity of dirt) spoon of detergent into tank.
C. Heat water with lid opened.
D. Pour out water after 30 minutes, and then wash the tank with clean water.
E. If cannot remove all dirt, please repeat process.
F. Residual dirt will not affect health, please feel free to use the machine.

6.4 Quarter Maintenance

For guaranteeing the purity and quality of distilled water, you need to replace the faucet regularly. The step as follows:
A. Push the left or right side of faucet, and then pull it, you can take it off.
B. According to the way of taking, you can put it on in an opposite way.
C. Repeat those steps when it needs to replace.

6.5 Power Off

Please turn off the power switch when no longer in use.
Please pull the end of power cord out of electric outlet when no longer in use.

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Chapter 1 Introduction

1.1 Attention

- ✦ This manual contains necessary and sufficient information to operate the system safely, like optimal use, a safe and reliable operation, regular and correct servicing requirements.
- ✦ Read and understand all instructions in this manual before attempting to use the product.
- ✦ Keep this manual with the equipment at all times. Periodically review the procedures for operation and safety precautions.

1.2 Application Range

- ✦ The medical distiller is designed by our company, which is for the places that needs High-purity water, such as clinic, Beauty Studio, hospital, laboratory.
- ✦ Operated by adult. Considering Ergonomic, it outstands on convenient using and safety, and continually, it can produce pure water 1L per hour in average.

1.3 Method

The pure water producing method is to transform liquid water into gaseous water by heating, then to achieve pure water (distilled water) through cooling.